-  **(EN) Enviro Swab**
-  **(FR) Enviro Swab**
-  **(DE) Umfeld-Tupfer**
-  **(IT) Tampone ambientale**
-  **(ES) Hisopo Enviro Swab**
-  **(NL) Enviro Swab**
-  **(SV) Enviro provtagningspinne**
-  **(DA) Enviro vatpind**
-  **(NO) Enviro Swab**
-  **(FI) Ympäristönäytteenottotikku**
-  **(PT) Enviro Swab**
-  **(EL) Στυλεός Enviro**
-  **(PL) Wacik do pobierania próbek Enviro**
-  **(HU) Enviroswab**
-  **(CS) Tampóny pro stěry z prostředí**
-  **(RO) Tampon Enviro**
-  **(RU) Enviro тампон**
-  **(TR) Enviro Pamuklu Çubuk**
-  **(JA) エンバイロスワブ 環境用拭取りスワブ**
-  **(ZH) Enviro 塗抹棒**

Product Instructions

Enviro Swab

PRODUCT DESCRIPTION AND INTENDED USE

A pre-moistened swab for environmental testing.

Neogen assumes that the Neogen® Enviro Swab be utilized by technicians that have been properly trained to follow microbiological techniques.

Neogen has not documented the Neogen Enviro Swab for use in industries other than food and beverage. For example, Neogen has not documented the Neogen Enviro Swab for testing water, pharmaceuticals or cosmetics. The Neogen Enviro Swab has not been evaluated with all possible surfaces or with all possible strains of bacteria.

SAFETY

The user should read, understand, and follow all safety information in the instructions for the Neogen Enviro Swab. Retain the safety instructions for future reference.

⚠ WARNING: Indicates a hazardous situation, which, if not avoided, could result in death or serious injury and/or property damage.

⚠ WARNING

Do not use the Neogen Enviro Swab in the diagnosis of conditions in humans or animals.

The user must train its personnel in current proper testing techniques: for example, Good Laboratory Practices or ISO/IEC 17025¹.

To reduce the risks associated with bacterial infection:

- Always follow standard lab safety practices, including wearing appropriate protective apparel while handling reagents and contaminated samples.
- Avoid contact with the contents of the Neogen Enviro Swab.
- Dispose of contaminated Neogen Enviro Swabs according to applicable Federal, State, and/or local government regulations and/or applicable laboratory procedures.

Consult the Safety Data Sheet for additional information.

For information on documentation of product performance, visit our website at www.neogen.com or contact your local Neogen representative or distributor.

USER RESPONSIBILITY

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.neogen.com, or contact your local Neogen representative or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results. The food sample itself may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any Neogen Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

LIMITATION OF WARRANTIES / LIMITED REMEDY

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, NEOGEN DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any Neogen Food Safety Product is defective, Neogen or its authorized distributor will, at its option, replace or refund the purchase price of the product. These are your exclusive remedies. You must promptly notify Neogen within sixty days of discovery of any suspected defects in a product and return it to Neogen. Please contact your Neogen representative or authorized Neogen distributor for any

further questions.

LIMITATION OF NEOGEN LIABILITY

NEOGEN WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall Neogen's liability under any legal theory exceed the purchase price of the product alleged to be defective.

STORAGE

When not in use, store Neogen Enviro Swabs between 2-8°C.

△ DISPOSAL

When testing is complete the Neogen Enviro Swab may contain microorganisms that may be a biohazard. Do not use the Neogen Enviro Swab past its expiration date. Expiration date and lot number are noted on the outside label of the box.


When testing is complete, follow current industry standards for the disposal of contaminated waste. Consult the Safety Data Sheet for additional information and local regulations for disposal.

INSTRUCTIONS FOR USE

Follow all instructions carefully. Failure to do so may lead to inaccurate results.

Use a fresh Enviro Swab for each sample area.

1. Label the Enviro Swab with sample identification.
2. Remove the tab from the top of the Enviro Swab (Figure 1).
3. Hold the tube firmly in one hand and release the swab by pushing the top of the tube with your thumb (Figure 2).
4. Gently push the tip of the swab to the side of the tube to remove excess diluent. Do not touch the swab or stem (Figure 3).
5. Carefully remove the swab from its tube and swab the area to be sampled (Figure 4).
6. Replace the swab in its tube and push firmly to lock the cap.

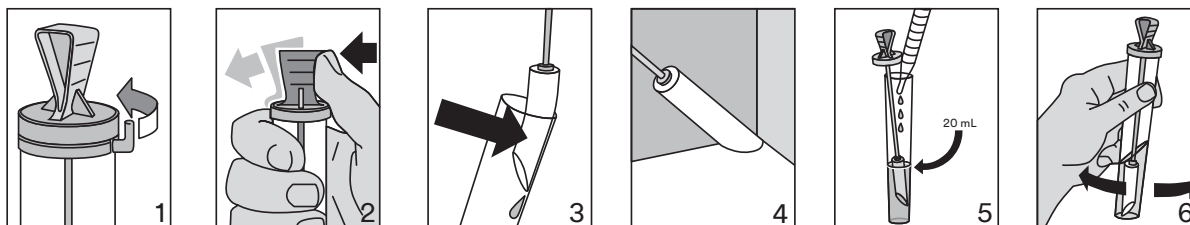7. Testing of the Neogen Enviro Swab must begin within 48 hours of sampling. Store the Neogen Enviro Swab at 2-8°C until testing begins.
8. Carefully lift the swab from its tube and aseptically pipette sufficient medium into the tube to cover the swab, approximately 20 mL (Figure 5).
9. Push the swab firmly back into the tube to lock the cap. Mix well by vortexing or shaking (Figure 6).
10. Incubate the Neogen Enviro Swab according to your method for environmental samples.

For further information refer to the Neogen® Enviro Swab "Interpretation Guide." If you have questions about specific applications or procedures, please visit our website at www.neogen.com or contact your local Neogen representative or distributor.

REFERENCES

1. ISO/IEC 17025. General requirements for the competence of testing and calibration laboratories.

Refer to the current versions of the standard methods listed above.

EXPLANATION OF SYMBOLS

info.neogen.com/symbols

Neogen Food Safety

Neogen Corporation

620 Leshar Place
Lansing, MI 48912 USA
neogen.com

Neogen Europe Ltd.

The Dairy School
Auchincruive
Ayr, KA6 5HU
Scotland, UK

Neogen Ireland, Ltd.

Bray Business Park, Bray
Co. Wicklow
A98YV29, Ireland

Neogen Corporation

620 Leshar Place Lansing, MI 48912 USA
www.neogen.com

Neogen is a trademark of Neogen Corporation.
© Neogen Corporation 2024. All rights reserved.
FS00878B

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