


INSTALLING DATA MANAGER SOFTWARE



To install Data Manager, run **setup.exe** from the installation CD. Be sure to check the **ENABLE RFID** box.

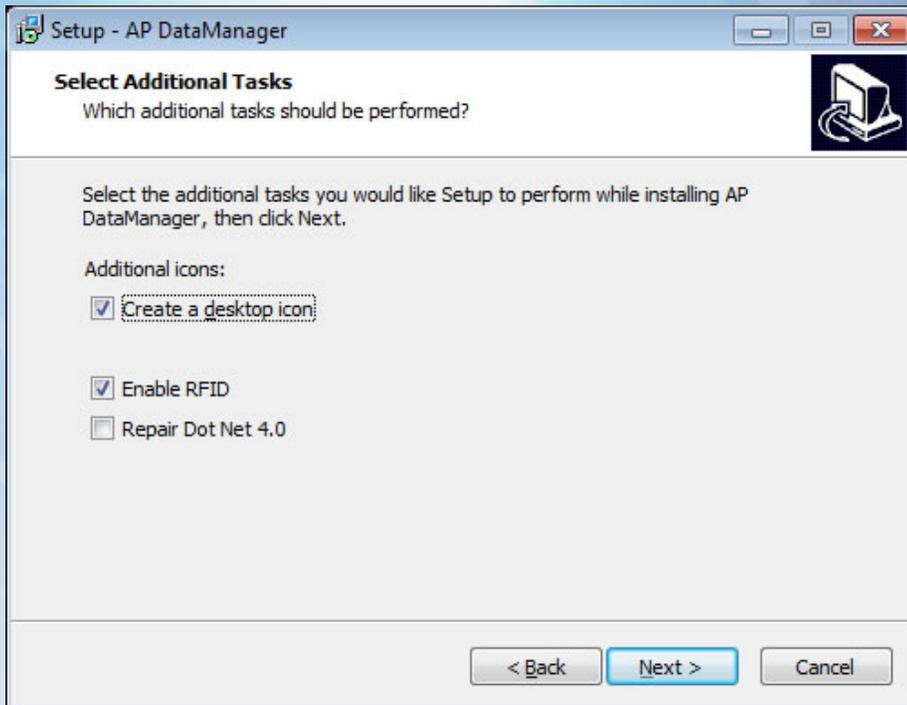

USA/Canada
800/234-5333 or 517/372-9200
foodsafety@neogen.com


Neogen Europe
+ 44 (0) 1292 525 640
0800 182 7721
info_de@neogeneurope.com

Neogen Latinoamerica
+ 52 (55) 5254-8235, + 52 (55) 5203-1198
informacion@neogenlac.com

Neogen do Brasil
+ 55 19 3935.3727
info@neogendobrasil.com.br

1 CREATE AND SYNC TEST PLAN

1. Use Data Manager to create a test plan for **Groups** and **Sites**.
 - a. **Groups** are organized on the left, **Sites** are organized on the right.
 - b. **Descriptions** can be added to each site.
 - c. Select the **Water** column to indicate a water sampler should be used during testing.
 - d. **Fail** threshold level can be set for each site.
 - e. **Marginal** threshold level can be set for each site.
 - f. **Mandatory site** can be identified.
 - g. **Select** or **Deselect** a site to be part of the synced plan.
 - h. Click on the **trash can** to delete a group or site.
2. Power on AccuPoint Advanced reader to the main screen and select **Sync to reader**.

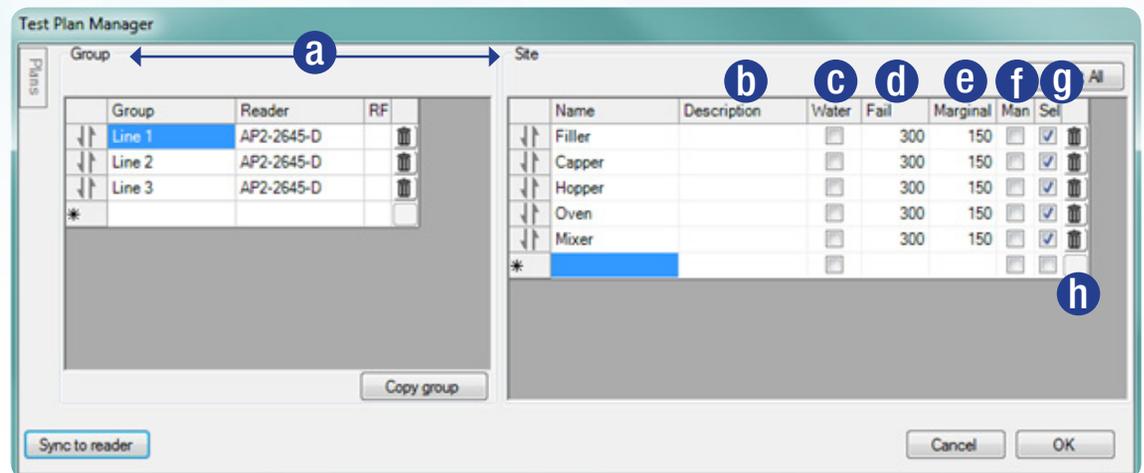

2 PROGRAM RFID TAGS



3. With the AccuPoint Advanced reader connected via USB, select **Program RFID Tags**.
4. Hold the antenna (on the back of the AccuPoint Advanced reader) near an RFID tag to assign the first group of sites.
5. Repeat the process for each group until all the tags have been assigned.
6. When all tags are assigned, select the  icon on the lower left of the display to accept the mapping.

3 UPLOAD RFID MAP



7. Reconnect AccuPoint Advanced reader to USB.
8. Select **Upload RFID Map**.

NOTE: An icon  will appear beside the Group when a RFID tag has been assigned. Use the **Clear RFID Map** command remove all RFID tags assigned to Groups.

PROGRAM MODE ICONS:

-  Select this to accept remapping of a RFID tag.
-  Select this to cancel the remapping of a RFID tag.
-  Indicates a group does not have a RFID tag assignment.
-  Indicates a group has a RFID tag assignment.
-  Shows when AccuPoint Advanced Reader is in RFID Program Mode.
-  Select to exit RFID Program Mode.



USA/Canada
800/234-5333 or 517/372-9200
foodsafety@neogen.com

Neogen Europe
+ 44 (0) 1292 525 640
0800 182 7721
info_de@neogeneurope.com

Neogen Latinoamerica
+ 52 (55) 5254-8235, + 52 (55) 5203-1198
informacion@neogenlac.com

Neogen do Brasil
+ 55 19 3935.3727
info@neogendobrasil.com.br