



T-1

PT-2

PT-3

Seek and Destroy *Listeria* Right Now.

Neogen® Molecular Detection Assay – *Listeria* Right Now™

Introducing a new test that delivers rapid results, no enrichment step required.

Optimize your workflow with faster, simpler testing by incorporating *Listeria* Right Now™ into your environmental monitoring program. Now, you can complete testing, read results, reclean, and retest all in the same shift, while receiving actionable results in about two hours from the standard 24-48 hours.

Rapid. Reliable. Right Now.

Now you have the ability to control contamination entry points in real time. Rapid results help manage entry points effectively and safeguard your production environment from *Listeria* species contamination.



Swab with Ease

Start with a sample collection device designed to efficiently swab surfaces up to 12"x12" and access narrow cracks and crevices.



Innovative Reagent Pellets

Each tube contains a blue, color-coded dehydrated pellet that reconstitutes easily when the sample is added and mixed by pipetting.



Results in ~2 Hours

No enrichment is required, allowing real-time decisions to be made during the same shift.

Neogen® Molecular Detection System

- User-friendly software with intuitive user interface for quick and easy setup
- Real-time results help make critical decisions faster
- Color-coded assays by organism and matching software screens allow for better tracking and easy identification
- Ability to run simultaneously with other qualitative assays
- Flexibility to test up to 96 samples in each run



Ordering Information

Description	Cat. No.	SKU	Contents
Neogen® Molecular Detection Assay – <i>Listeria</i> Right Now™	MDALRN96	700007660	96 Tests/Kit
Neogen® Molecular Detection Right Now™ Matrix Control	RNMC96	700007661	96 Tests/Kit
Neogen® Molecular Detection Instrument and Accessory Kit	MDS100	700002195	1 Each
Neogen® Molecular Detection Heat/Chill Block Insert	MDSHBIN	700000145	1 Each

Learn more at: info.neogen.com/ListeriaRightNow



Neogen Corporation, 620 Leshar Place, Lansing, MI 48912 USA.

© Neogen Corporation 2025. All rights reserved. Neogen is a registered trademark of Neogen Corporation.

All other marks mentioned are the property of their respective owners.

FS02116_0325