-  (EN) Quick Swab
-  (FR) Quick Swab
-  (DE) Quick Swab
-  (IT) Tampone rapido
-  (ES) Quick Swab
-  (NL) Quick Swab
-  (SV) Quick Swab
-  (DA) Quick Swab
-  (NO) Quick Swab
-  (FI) Quick Swab Näytetikku
-  (PT) Quick Swab
-  (EL) Στυλεός Δειγματοληψίας Γρήγορης Λήψης
-  (PL) Szybka wymazówka
-  (HU) Quick Swab gyors kenetvevő
-  (CS) Rychlý stěr
-  (RO) Tampon de recoltare rapidă
-  (RU) Тампон для отбора образцов «Quick swab»
-  (TR) Hızlı Svab
-  (JA) クイック・スワブ
-  (ZH) 快速涂抹棒
-  (TH) ควิกสวอป
-  (KO) 퀵 스왑

Product Instructions

Quick Swab

Product Description and Intended Use

The Neogen® Quick Swab is a ready-to-use environmental swab system that can be used wet or dry, and it is intended to be used in the food and beverage industry for surface sampling. Neogen Quick Swab is designed to be used in conjunction with any Neogen® Petrifilm® Plate or other culture media. The Neogen Quick Swab uses a sodium bisulfite-free letheen broth to facilitate the recovery of bacteria and will deliver approximately 1.0 mL of sample if used in the prescribed manner.

Neogen Quick Swabs are aseptically filled, and gamma irradiated.

Safety

The user should read, understand, and follow all safety information in the instructions for the Neogen Quick Swab. Retain the safety instructions for future reference.

⚠ CAUTION Indicates a hazardous situation, which, if not avoided, could result in minor or moderate injury and/or property damage.

⚠ CAUTIONS

- The Neogen Quick Swab does not contain any indicator dyes. Therefore, it alone will not detect bacterial colonies.
- The Neogen Quick Swab must be used with a plating media such as Neogen Petrifilm Plates.
- The Neogen Quick Swab has not been tested with all possible food processing sanitation products.
- Do not use the Neogen Quick Swab in the diagnosis of conditions in humans or animals.
- Do not use the Neogen Quick Swab more than once.
- The user must train its personnel in proper testing techniques, for example: Good Laboratory Practices⁽¹⁾ ISO 7218⁽²⁾ or ISO 17025⁽³⁾.
- Consult the Safety Data Sheet for additional information.
- For information on documentation of product performance, visit our website at www.neogen.com or contact your local Neogen representative or distributor.

User Responsibility

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.neogen.com, or contact your local Neogen representative or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results. The food sample itself may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any Neogen Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

Limitation of Warranties / Limited Remedy

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, NEOGEN DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any Neogen Food Safety Product is defective, Neogen or its authorized distributor will, at its option, replace or refund the purchase price of the product. These are your exclusive remedies. Please contact your Neogen representative or authorized Neogen distributor for any further questions.

Limitation of Neogen Liability

NEOGEN WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall Neogen's liability

under any legal theory exceed the purchase price of the product alleged to be defective.

Storage

Store Neogen Quick Swab at temperatures $\leq 8^{\circ}\text{C}$ ($\leq 46^{\circ}\text{F}$). Exposing the Neogen Quick Swab to temperatures $> 25^{\circ}\text{C}$ ($> 77^{\circ}\text{F}$) can affect the performance of the product. Lot numbers and expiration dates are noted on each box or case of Neogen Quick Swabs as well as on each swab.

Δ Disposal

As with all environmental surface sampling swabs, once used, the Neogen Quick Swab may be a potential biohazard. Follow current industry standards for disposal.


Instructions for Use

Follow all instructions carefully. Failure to do so may lead to inaccurate results.


Prior to use, visually inspect each Neogen Quick Swab. Use only if the letheen broth is in the bulb end of the swab and the color of the broth is clear amber, not turbid or cloudy.

Figure 1


1 mL Plating Instructions - Wet Swabbing Method -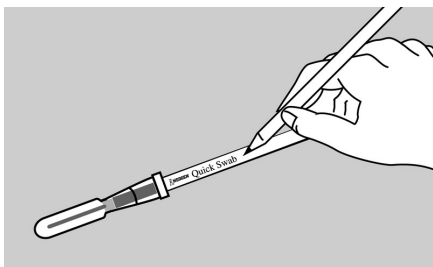


1 Label each Neogen Quick Swab.

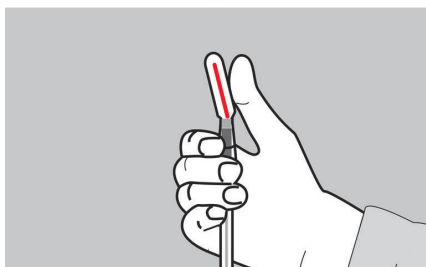
2 Bend the red snap valve at a 45° angle until you hear the valve break. This allows the letheen broth to flow into the tube end and wet the swab bulb.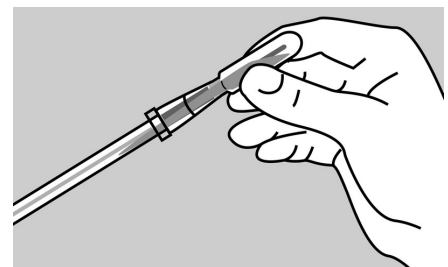


3 Squeeze the bulb of the swab to transfer all of the letheen broth to the tube end of the swab.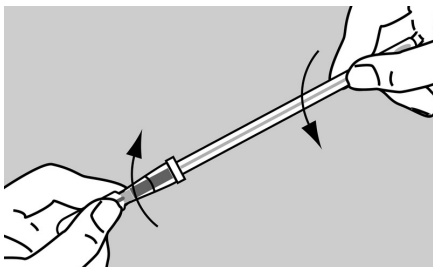




4 Twist and pull apart the bulb end of the swab from the tube end of the swab that contains the letheen broth.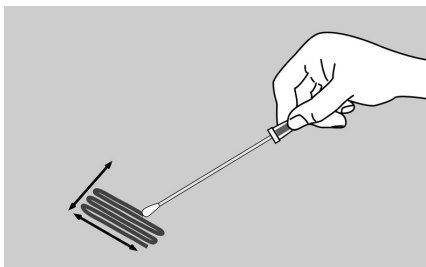

5 Hold the swab handle to make a 30° angle with the surface. Rub the swab slowly and thoroughly over the desired surface area. Rub the swab three times over this surface, reversing direction between alternating strokes.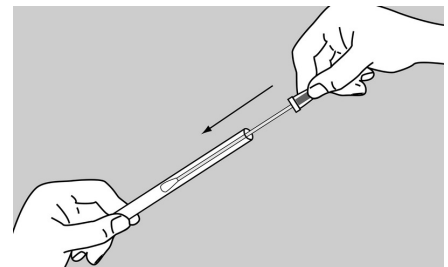

6 After sampling is complete, securely insert the swab back into the swab tube and transport to the lab for inoculation. Plate the letheen broth swab solution as soon as possible.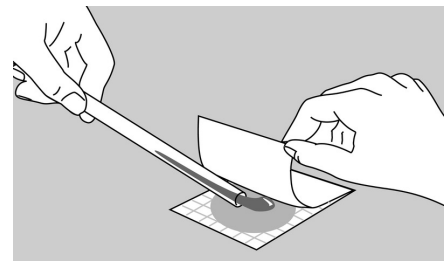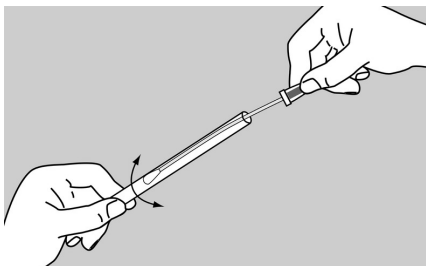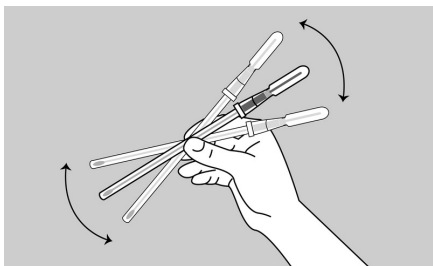


7 In the lab, vigorously shake or vortex the Neogen Quick Swab for 10 seconds, to release bacteria from the swab tip.

8 Wring out the contents of the swab tip by pressing and twisting the swab against the wall of the tube.


9 Carefully pour the entire contents of the tube onto a Neogen Petrifilm Plate. Follow current industry standards for disposal.

Figure 2


1 mL Plating Instructions - Dry Swabbing Method -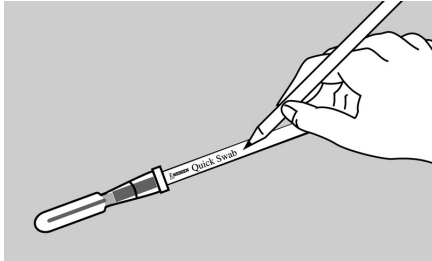


1 Label each Neogen Quick Swab.

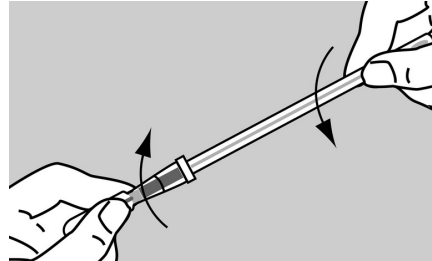
2 Twist and pull apart the bulb end of the swab from the tube end of the swab that contains the letheen broth.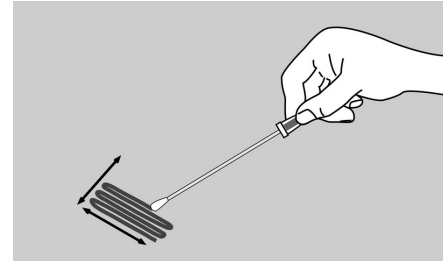


3 Hold the swab handle to make a 30° angle with the surface. Rub the swab slowly and thoroughly over the desired surface area. Rub the swab three times over this surface, reversing direction between alternating strokes.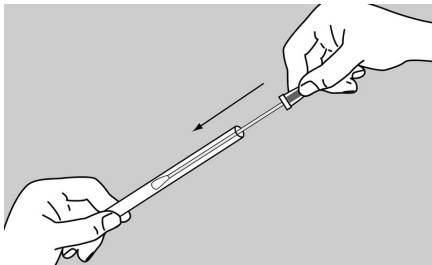

4 After sampling is complete, securely insert the swab back into the swab tube and transport to the lab for inoculation. Plate the letheen broth swab solution as soon as possible.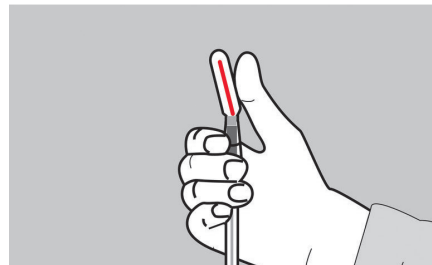

5 Bend the red snap valve at a 45° angle until you hear the valve break. This allows the letheen broth to flow into the tube end and wet the swab.

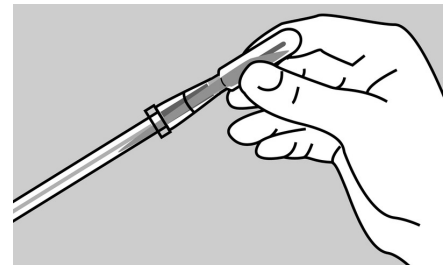6 Squeeze the bulb of the swab to transfer all of the letheen broth to the tube end of the swab.

To complete inoculation, continue with steps 7-9 (See Figure 1 above)

Multiple mL Plating Instructions:

Plating multiple bacteria tests using the Neogen Quick Swab (Wet or Dry Method)

- a. Complete Steps 1 through 6 (Figures 1 or 2 above) for either the Wet or Dry swabbing method.
- b. Remove the Neogen Quick Swab from the tube. Add 1-3 mL of sterile diluent to swab tube. Replace the swab in the tube.
- c. To prepare the sample for inoculation, vigorously shake or vortex Neogen Quick Swab for 10 seconds to release the bacteria from the swab tip.
- d. Wring out the contents of the swab tip by pressing and twisting the swab against the wall of the tube.
- e. Use your thumb to bend the swab tube at a 90° angle at the highest mark that has a diluent above it. Pour off a 1 mL aliquot onto a Neogen Petrifilm Plate. Repeat until the entire sample is used.
- f. Follow current industry standards for disposal.

The entire contents of Neogen Quick Swab may also be added to any sterile dilution blank. Pipette 1 mL aliquots from the dilution blank onto Neogen Petrifilm Plates.

If used with Neogen Petrifilm Plates:

Refer to the appropriate Neogen Petrifilm Plate “Product Instructions.”

If you have questions about specific applications or procedures, please visit our website at www.neogen.com or contact your local Neogen representative or distributor.

NOTE: As with any swab sampling method, the best microbial practice is to re-sanitize the swabbed area after the swab sample has been taken.

References

1. U.S. Food and Drug Administration. Code of Federal Regulations, Title 21, Part 58. Good Laboratory Practice for Nonclinical Laboratory Studies.
2. ISO 7218. Microbiology of food and animal feeding stuffs – General requirements and guidance for microbiological examinations.
3. ISO/IEC 17025. General requirements for the competence of testing and calibration laboratories.

Refer to the current versions of the standard methods listed above.

Explanation of Symbols

info.neogen.com/symbols

Neogen Food Safety

Neogen Corporation
620 Leshar Place
Lansing, MI 48912 USA
neogen.com

Neogen Europe Ltd.
The Dairy School
Auchincruive
Ayr, KA6 5HU
Scotland, UK

Neogen Ireland, Ltd.
Bray Business Park, Bray
Co. Wicklow
A98YV29, Ireland

Neogen Corporation
620 Leshar Place Lansing, MI 48912 USA
www.neogen.com

Neogen and Petrifilm are trademarks of Neogen Corporation.
© Neogen Corporation 2024.
All rights reserved.
FS00831B

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