Read instructions carefully before using

Allergen Environmental Swabbing Kit

INTENDED USE

The 100 swabs and dropper tips in this kit are intended to be used to sample environmental surfaces for the possible presence of food allergen residues. The swabs allow for the sampling of all potential food contact surfaces, including all types of processing and storage equipment, as well as conveyors and work surfaces. The extraction solution required for use with the swabs for testing dry surfaces is provided with Neogen's Alert® and Reveal® food allergen screening test kits. Samples obtained from the use of this kit are intended to be analyzed with Neogen's screening test kits.

STORAGE

This kit may be stored at room temperature (18–30°C, 64–86°F).

STORAGE OF ENVIRONMENTAL SWAB SAMPLES

If environmental swabs are to be stored or shipped, they should be moistened with 100–200 µL of extraction buffer at 4–8°C and also should be shipped under these conditions to preserve the samples' integrity. Due to diminished protein solubility over time, Neogen recommends testing within 24 hours from the time of collection. This is considered optimum for peak recovery of allergenic residues. In situations where this is not logistically feasible, the samples can be analyzed up to 72 hours from the time of collection with valid results.

MATERIALS PROVIDED

1. 100 sterile swabs
2. 100 disposable dropper tips

MATERIALS RECOMMENDED BUT NOT PROVIDED


1. Bottle to prepare extraction solution, 1 L (Neogen Item 9472)
2. Bottle to prepare working solution, 150–250 mL
3. Extraction solution appropriate for specific food allergen test kit
4. Ethanol, 190 proof (for gliadin testing only)
5. Distilled or deionized water
6. Vortex mixer, or equivalent (Neogen item 9494)
7. Food allergen test kit

SAMPLE PREPARATION AND EXTRACTION

For peanut, almond, crustacea*, hazelnut, lupine, egg, milk, mustard, soy or soy flour

Test samples should be collected according to accepted sampling techniques (if necessary, see Neogen's Food Allergen Handbook). The extraction solution needed for this procedure varies with the target food allergen, and is supplied with the Alert or Reveal food allergen test kit. Environmental swabs should be extracted and analyzed within **4 hours** of collection. See **Storage of environmental swab samples** for extended storage and shipping instructions.

1. Follow the instructions in an Alert or Reveal test kit to prepare a target allergen specific extraction solution.
2. Prepare a working solution by combining 125 mL of the prepared extraction solution in a separate container, and add 1 level scoop of extraction additive from the test kit to the container. Prepare a fresh working solution daily. **NOTE:** Prepare 125 mL of working solution for every 25 environmental swabs to be tested.
3. Preheat the working solution to 60°C (140°F)*. Shake container to dissolve additive.
4. Gather the sample with a swab, using one of the following methods:
For dry surfaces: Open a new swab and wet with extraction solution. Swab a 10 x 10 cm area by using a crosshatch technique. **NOTE:** Do not use working solution to moisten swabs.
For wet surfaces: Open a new swab and swab a 10 x 10 cm area by using a crosshatch technique. Do not moisten swab prior to use.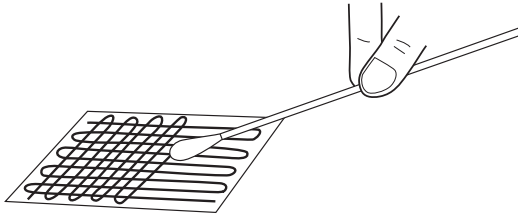

5. Return the swab to its original tube once sampling is complete. Remember to label each tube.
6. Remove the swab from its tube, and add 5 mL of 60°C* working solution to the tube. Mix by placing the swab back into the tube and shaking for **2 minutes** by hand (inverting tube), or for **30 seconds** with a vortex mixer.
7. Remove the swab from its tube.
8. **Alert:** If using an Alert test kit, place a new sample dropper tip onto the tube. The solution in the tube now serves as the sample. Continue sample analysis by following testing instructions in the Alert allergen test kit. **NOTE:** Use caution when inverting the tube, as some liquid may drip.
Reveal: If using a Reveal test kit, transfer the appropriate volume of the supernatant to a sample tube. Continue sample analysis by following testing instructions in the Reveal allergen test kit.

***Crustacea allergen swabs should be extracted in 30°C (86°F) working solution.**

SAMPLE PREPARATION AND EXTRACTION

For gliadin (non-heat processed samples)

NOTE: Contact Neogen Technical Services for the gliadin cocktail procedure for heat processed samples.

Test samples should be collected according to accepted sampling techniques (if necessary, see Neogen's Food Allergen Handbook). Environmental swabs should be extracted and analyzed within **4 hours** of collection. See **Storage of environmental swab samples** for extended storage and shipping instructions.

For gliadin R5 kits (Neogen items 8510 and 8511): These kits require the use of a 60% ethanol solution. To prepare, combine 6 parts of 190 proof ethanol with 4 parts distilled water.

For gliadin kits (Neogen items 8480 and 8481): These kits require the use of a 40% ethanol solution. To prepare, combine 4 parts of 190 proof ethanol with 6 parts distilled water.


1. Prepare appropriate ethanol solution (see above).
2. Prepare a working solution by placing 20 mL of the appropriate ethanol solution in a clean container. Add 1 scoop extraction additive from the test kit and swirl to mix.

NOTE: Prepare 20 mL of working solution for every four environmental swabs to be tested.

3. Gather sample with a swab, using one of the following methods:

For dry surfaces: Open a new swab and wet with the appropriate ethanol solution. Swab a 10 x 10 cm area by using a crosshatch technique. **NOTE:** Do not use working solution to moisten swabs.

For wet surfaces: Open a new swab and swab a 10 x 10 cm area by using a crosshatch technique. Do not moisten swab prior to use.

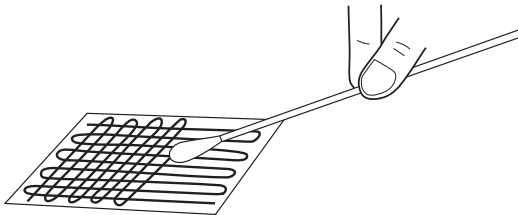4. Return the swab to its original tube once sampling is complete. Remember to label each tube.
5. Remove the swab from its tube, and add 5 mL of ethanol solution with additive to the tube. Mix by placing the swab back into the tube and shaking for **2 minutes** by hand (inverting tube), or for **30 seconds** with a vortex mixer.
6. Remove the swab from its tube and place a new dropper tip onto the tube.
7. Dilute by adding 3 drops of extract into a prefilled sample dropper bottle. (Provided with the appropriate test kit.)

CUSTOMER SERVICE

Neogen Customer Assistance and Technical Services can be reached by using the contact information on the back of this booklet. Training on this product, and all Neogen test kits, is available.

MSDS INFORMATION AVAILABLE

Material safety data sheets (MSDS) are available for this test kit, and all of Neogen's test kits, on Neogen's website at www.neogen.com, or by calling Neogen at 800/234-5333 or 517/372-9200.

TERMS AND CONDITIONS

For Neogen's full terms and conditions, please visit www.neogen.com/Corporate/termsconditions.html.

WARRANTY

Neogen Corporation makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, Neogen will provide a replacement of the product. Buyer assumes all risk and liability resulting from the use of this product. There is no warranty of merchantability of this product or of the fitness of the product for any purpose. Neogen shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

TEST KITS AVAILABLE FROM NEOGEN

Natural toxins

- Aflatoxin, DON, ochratoxin, zearalenone, T-2/HT-2 toxins, fumonisin, histamine

Foodborne bacteria

- *E. coli* O157:H7, *Salmonella*, *Listeria*, *Listeria monocytogenes*, *Campylobacter*, *Staphylococcus aureus*

Sanitation

- ATP, yeast and mold, total plate count, generic *E. coli* and total coliforms, protein residues

Food allergens

- Almonds, crustaceans, eggs, gliadin, hazelnut, lupine, milk, mustard, peanuts, sesame, soy, walnuts

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