Petrifilm®

Important Information

- ① **Automated Feeder**
- ② **Chargeur automatisé**
- ③ **Automatisierter Verteiler**
- ④ **Caricatore automatico**
- ⑤ **Alimentador automatizado**
- ⑥ **Geautomatiseerde feeder**
- ⑦ **Automatisk matare**
- ⑧ **Automatiseret forsyner**
- ⑨ **Automatisert forsyner**
- ⑩ **Syöttöautomaatti**
- ⑪ **Alimentador automatizado**
- ⑫ **Αυτοματοποιημένο σύστημα τροφοδοσίας**
- ⑬ **Automatyczny podajnik**
- ⑭ **Устройство автоматической подачи**
- ⑮ **Otomatik Besleyici**
- ⑯ **自動フィーダー**
- ⑰ **自动送料器**
- ⑱ **เครื่องป้อนอัตโนมัติ**
- ⑲ **자동 공급 시스템**

Petrifilm®

Important Information

Intended Use

The Neogen Petrifilm® Automated Feeder processes Neogen Petrifilm® Plates and is not intended to be used with any other manufactured products. The Neogen Petrifilm Automated Feeder utilizes vacuum and mechanical forces to transfer up to 300 Neogen Petrifilm Plates from an input bin into the Neogen Petrifilm Plate Reader Advanced. Once processed by the Neogen Petrifilm Plate Reader Advanced, the plates will be transferred to a sorting station in which the Neogen Petrifilm Automated Feeder will transfer each plate to one of two output bins based on user defined criteria. The Neogen Petrifilm® Plate Manager software is designed for the transfer, organization, data storage and data management for the results of the Neogen Petrifilm Plates.

Neogen has not designed nor documented the Neogen Petrifilm Automated Feeder, and its power supply, or Neogen Petrifilm Plate Manager for use with other manufacturers' products. Use with other manufacturers' designs or use in any other application that has not been evaluated or documented by Neogen may lead to an unsafe condition. Although the Neogen Petrifilm Automated Feeder system is accurate, all results should be reviewed and confirmed manually by the user. The Neogen Petrifilm Automated Feeder system is not intended to be used in the diagnosis of conditions in humans or animals. The user is responsible to know and follow applicable workplace regulations.

Neogen anticipates that the Neogen Petrifilm Plates, equipment and software will be utilized by technicians that have been properly trained on the Neogen Petrifilm Plate test methods, the Neogen Petrifilm Automated Feeder, Neogen Petrifilm Plate Reader Advanced and the Neogen Petrifilm Plate Manager.

If this product is used in a manner that is not specified, the protection provided by the product may be impaired.

Neogen is certified to ISO (International Organization for Standardization) 9001 for design and manufacturing.

Safety Information

Please read, understand, and follow all safety information contained in these instructions prior to use of this Neogen Petrifilm Automated Feeder. Retain these instructions for future reference.

User Responsibility

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.neogen.com, or contact your local Neogen representatives or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any Neogen product do not constitute a guarantee of the quality of the matrices or processes tested.

Explanation of Safety Related Symbols	
	Warning: Hazardous Voltage
	Attention: Read Accompanying Documentation
	Caution: Consult the instructions for use
	This product contains recyclable parts as per Article 219(5) of Italian Legislative Decree No. 152 of 2006. For information on recycling - please contact your nearest Neogen Service Center for advice.
	WEEE mark - This product contains electrical and electronic components and must not be disposed of using standard refuse collection. Please consult local directives for disposal of electrical and electronic equipment.
	TUV NRTL mark is a conformity marking that indicates conformity with the applicable electrical safety and EMC/EMI requirements for products sold within the U.S. and Canada.
	UK Conformity Assessed (UKCA) marking is a conformity mark that indicates conformity with all applicable electrical product safety and EMC/EMI requirements for products sold within Great Britain
	Japan's Ministry of Economy, Trade and Industry (METI) established the Electrical Appliance and Material Safety Law (DENAN) to regulate the safety of electronic and electrical products.
	National Institute of Metrology, Standardization and Industrial Quality - Inmetro governs the (Inmetro) marking as a conformity mark that indicates conformity with the applicable safety requirements for products sold within Brazil.
	The RoHS mark indicates products sold within the European Economic Area are compliant to RoHS Directive 2011/65/EU. Therefore this electrical and electronic equipment (EEE) dose not contain restricted substances cited in the RoHS 3 directive.
	The Korean Ministry of Trade, Industry and Energy (MOTIE) governs the Korea Certification Mark (KC Mark) as a conformity mark that indicates compliance with the applicable electrical safety requirements, EMC, and energy efficiency for products sold within S. Korea.
	CE mark (conforms to applicable requirements for European Economic Area).
	RCM mark (Australia and New Zealand electrical safety and EMC)
	Date of manufacture YYYY-MM-DD

Explanation of Signal Word Consequences

- ⚠ WARNING:** Indicates a hazardous situation which, if not avoided, could result in death or serious injury.
- ⚠ CAUTION:** Indicates a hazardous situation which, if not avoided, could result in minor or moderate injury and/or property damage.

⚠ WARNING

To reduce the risk associated with hazardous voltage:

- Use indoors only.
- Do not spill liquids on device or immerse in liquids.
- The Personal Computer (PC) that the Neogen Petrifilm Automated Feeder connects to must be an approved certified device (mark with UL, CSA, ETL, TUV, or other required regulatory safety agency in your region) and meet requirements specified in the user manual.
- The USB cable must only be connected to an approved listed/certified device.

To reduce the risk of serious injury

- In order to avoid injury the Petrifilm Automated Feeder should not be lifted with out the use of a forklift or lift table.

⚠ CAUTION

To reduce the risk associated with EMC interference:

- Modifications to this device or power supply shall not be made without the written consent of Neogen. Unauthorized modifications may void the authority granted under federal communication rules permitting the operation of this device.

To reduce the risk associated with misinterpretation of results:

- Follow all instructions provided in the Neogen Petrifilm interpretation guides, Petrifilm product instructions and the Neogen Petrifilm Plate Reader Advanced User Manual and the Automated Feeder User Manual.
- In order to avoid inaccurate results users should only use the Petrifilm Automated Feeder with the Petrifilm Plate tests that are listed in the Neogen Petrifilm Plate Reader Advanced user manual.

To reduce risk associated with bacterial infection and workplace contamination:

- In order to avoid exposure to microorganisms, the user should clean the instrument in the event damage is done to the Petrifilm Plate resulting in media being exposed to the environment.
- In order to avoid exposure to excessive dust and powders the user should perform routine cleaning of the instrument.
- In order to avoid exposure to microorganisms the user should follow good laboratory practices.
- In order to avoid cross contamination users should follow good laboratory practices.

To reduce the risk of injury, instrument damage or damage to the Neogen Petrifilm Plates:

- In order to avoid injury user should not operate the Petrifilm Automated Feeder when the bin doors are open.

- In order to avoid injury, the user should notify service personnel if the primary door is not operating as expected.
- In order to avoid the Petrifilm Automated Feeder jamming, the Petrifilm Plates should be placed in the bin so that the hinged side of the Petrifilm Plate feeds into the PPRA first.

To reduce the risks associated with exposure to biohazards and environmental contamination:

- Local laws and regulations should be followed for the disposal of hazardous materials.

Electrically Safe Work Condition

- Exposed energized equipment and parts will first be deenergized, locked/tagged out, tested to verify that there is no voltage, and grounded if necessary to protect workers, before any maintenance or repair work is performed. Only a qualified person will deenergize, lock/tag out, and test electrical parts and equipment.

Protective Earthing

- Installed power cord assembly must have a protective-earth conductor. Protective earthing connections shall be such that disconnection of a protective earth at one point in a unit or a system does not break the protective-earthing connection to other parts or units in a system, unless the relevant hazard is removed at the same time.

General Safe Work Practices

- All employees working on or near electrical equipment will follow general safe work practices, including:
 - Maintain good housekeeping procedures.
 - Plan and analyze safety in each step of a project.
 - Document work.
 - Use properly rated test equipment and verify its condition and operation before and after use.
 - Practice applicable emergency procedures.
 - Maintain electrical equipment in accordance with the manufacturer's instructions.
 - Plan work projects through an approved work control process.

Grounding-type Equipment

- A flexible cord used with grounding-type equipment will contain an equipment grounding conductor. Attachment plugs and receptacles may not be connected or altered in a manner that would prevent proper continuity of the equipment grounding conductor at the point where plugs are attached to receptacles. Additionally, these devices may not be altered to allow the grounding pole of a plug to be inserted into slots intended for connection to the current-carrying conductors. Adapters which interrupt the continuity of the equipment grounding connection may not be used.

Electrical Maintenance and Repair Operations

- Only qualified personnel will perform repair or maintenance work on electrical conductors or circuits. If an electrical hazard is discovered while repairs or maintenance work is performed, any further work must be suspended until the hazard is addressed and corrective actions instituted.
- Qualified personnel performing such tasks as electrical repairs, modifications, and tests on energized conductors and circuits, parts, and equipment will comply with the following work practices.

Utilization Equipment

- Utilization equipment is subject to the same approval and acceptance requirements as that of electrical equipment. To be acceptable for installation and use, utilization equipment will be listed or labeled by a nationally recognized testing laboratory. Utilization equipment that is not listed or labeled will meet one of the requirements of 29 CFR 1910.399, Acceptable, (i) (ii), or (iii). Utilization equipment that is not listed or labeled will be examined, accepted, and documented by a qualified person. Utilization equipment will be used in accordance with its listing and labeling requirements.

Regulatory Compliance

HEALTH AND ELECTRICAL SAFETY INFORMATION

The Neogen Petrifilm Automated Feeder complies with the following standards as demonstrated by the CB Scheme Certificate and test report issued by TÜV SÜD America Inc. (UL): • UL/IEC/EN 61010-1:2010. Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements • Neogen Petrifilm Automated Feeder is TUV listed and carries the TUV mark with adjacent indicators “C” and “US” based on compliance to the standards UL 61010-1, CAN/CSA 22.2 No. 61010-1, (USA and Canada). The Neogen Petrifilm Automated Feeder complies with the CE mark related to the Low Voltage Directive (LVD) 2014/35/EU as confirmed in the Declaration of Conformity. The Neogen Petrifilm Automated Feeder complies with the RoHS Directive, Directive 2023/143 of the European Parliament and of the Council of 4 May 2023 on the restriction of the use of certain hazardous substances in electrical and electronic equipment. Neogen Petrifilm Automated Feeder complies with China RoHS II requirements. Neogen Petrifilm Automated Feeder complies with the WEEE Directive, Directive 2012/19/EU of the European Parliament and of the Council of 04 July 2012 on waste electrical and electronic equipment (WEEE).

Circuit protection is supported by the following primary safety critical component:

Component	Mfr Name	Model/ Type	Voltage (VAC)	Current (A)	Safety Listing
Fuse	Littelfuse	218 Series	250	10	UL 62368-1

EMC Compliance

The Neogen Petrifilm Automated Feeder complies with the following EMC standards as confirmed in the Certificate of Compliance generated by Neogen:

- IEC 61326-1:2012/EN 61326:2013 Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements
- EMC requirements of the CE mark EMC Directive 2014/30/EU. This equipment is not intended for use in residential environments and may not provide adequate protection to radio reception in such environments. The Neogen Petrifilm Automated Feeder complies with the Australian and New Zealand electrical safety and electromagnetic compatibility requirements as confirmed in the Supplier’s Declaration of Conformity that is linked to the Australian/ New Zealand Regulatory Compliance Mark (RCM). Innovation, Science and Economic Development Canada ICES-003 Compliance Label: CAN ICES-3 (A)/NMB-3(A). This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation. Note: This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial

environment. This equipment generates, uses, and can radiate radiofrequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at their own expense.

For information on documentation of product performance, visit our website at www.neogen.com or contact your local Neogen representative or distributor.

ENVIRONMENTAL OPERATING CONDITIONS

Environmental Condition	Operating Condition	Units
Indoor Use only	- only connect the data connections to a computer that conforms to IEC/EN/UL/CSA 60950-1 or 62368-1	
Altitude	2500 maximum	meters
Operating Temperature Range	15 to 35	°C
Relative Humidity	10 – 85 noncondensing	%
Storage Temperature Range	-20 to +60 (-4 to +140)	°C (°F)
Overvoltage	Category II	

Model # 700006589

AC Voltage for power supply Mains Supply voltage fluctuations up to ± 10% of nominal voltage Temporary overvoltages occurring on the mains supply	100-120	Volts
Frequency	50/60	Hertz
AC Current for power supply	8	Amps

Model # 700006590

AC Voltage for power supply Mains Supply voltage fluctuations up to ± 10% of nominal voltage Temporary overvoltages occurring on the mains supply	220-240	Volts
Frequency	50/60	Hertz
AC Current for power supply	4	Amps

INSTRUMENT SPECIFICATIONS

Dimensions	Specification	Units
Length	1182	mm
Width	486	mm
Height	781	mm
Weight	110	Kg
External Connectors	USB 2.0 Type B and 4 Pin Lock DIN	

FACTORY/MANUFACTURER

Factory Name	Levison Enterprises
Address	4470 Moline Martin Road
Locale	Millbury, OH 43447
Phone #	(419) 838-7365
Website Address	https://levisonenterprises.com/

Cleaning and Decontaminating the Neogen Petrifilm Automated Feeder

Lightly spray a disposable towel with 70% Isopropyl alcohol (IPA) or 70% Ethanol and gently wipe the outer and inner surfaces of the instrument. The disposable towel should be modestly moist and not dripping. Avoid the power and USB cable connection while cleaning the exterior surfaces.

Note: Please refer to the Cleaning and Maintenance section of the software for video tutorials on specific components to clean.

Limited Warranty

Neogen Hardware (“Hardware”) includes the Neogen Petrifilm Automated Feeder and any associated Software and other components supplied by Neogen which are described in the applicable Installation and User Manuals (“User Documentation”).

Limited Warranty and Disclaimer

Neogen warrants that for one year from the date of shipment that the Hardware will substantially perform in accordance with the User Documentation. THIS WARRANTY IS VOID IF: (A) THE HARDWARE HAS BEEN REPAIRED BY PERSONS NOT AUTHORIZED BY NEOGEN; OR (B) THE HARDWARE HAS BEEN ALTERED, MODIFIED, OR MISUSED; OR (C) THE HARDWARE IS USED WITH PRODUCTS, SUPPLIES, COMPONENTS OR SOFTWARE NOT SUPPLIED BY NEOGEN FOR USE WITH THE HARDWARE; OR (D) THE HARDWARE OR A COMPONENT IS USED FOR OTHER USES (FOR EXAMPLE USE WITH OTHER CIRCUIT BOARDS OR SOFTWARE) OR (E) THE HARDWARE HAS NOT BEEN MAINTAINED OR USED IN ACCORDANCE WITH THE USER DOCUMENTATION. UNLESS PROHIBITED BY LAW, THIS WARRANTY IS MADE IN LIEU OF ALL OTHER WARRANTIES, EXPRESS OR IMPLIED, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTY OF FITNESS FOR A PARTICULAR PURPOSE, THE IMPLIED WARRANTY OF MERCHANTABILITY, OR ANY IMPLIED WARRANTY ARISING OUT OF A COURSE OF DEALING OR OF PERFORMANCE, CUSTOM OR USAGE OF TRADE. NEOGEN DOES NOT WARRANT THAT THE HARDWARE AND SOFTWARE WILL FUNCTION ERROR FREE. If within one year after shipment, the Hardware does not conform to the express warranty set forth above, Neogen’s sole obligation and User’s sole remedy shall be, at Neogen’s option: 1) to repair or replace the non-conforming component; or, 2) refund the purchase price.

Limitation of Liability

UNLESS PROHIBITED BY LAW, NEOGEN WILL NOT BE LIABLE TO USER OR OTHERS FOR ANY OTHER DIRECT, INDIRECT, CONSEQUENTIAL, INCIDENTAL OR SPECIAL DAMAGES INCLUDING, FOR EXAMPLE, LOST PROFITS, BUSINESS, INVESTMENTS, OR OPPORTUNITIES EVEN IF NEOGEN HAS BEEN ADVISED OF THE POSSIBILITY OF SUCH DAMAGES. The parties agree that Neogen’s total cumulative liability to User for direct damages for all causes shall not exceed One Hundred Dollars, or the price paid for the Hardware, whichever is higher. Some states or countries may have laws which require liability rights different from those stated above. In such state or countries, the minimum required liability terms shall apply.

Disclaimer: Neogen disclaims all express and implied warranties including warranties of merchantability or fitness for a particular use. If product is defective, the exclusive remedy is, at Neogen's option, replacement, repair or refund. Except where prohibited by law, Neogen will not be liable for any further loss or damage arising from use of this product.

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.neogen.com.

Neogen Food Safety

Neogen Corporation

620 Lesher Place
Lansing, MI 48912 USA
Neogen.com

Neogen Europe Ltd.

The Dairy School
Auchincruive
Ayr, KA6 5HU
Scotland, UK

Neogen Ireland, Ltd.

Bray Business Park, Bray
Co. Wicklow
A98YV29, Ireland

Neogen Corporation

620 Lesher Place
Lansing, MI 48912 USA
www.neogen.com

Neogen is a trademark of Neogen Corporation.
© Neogen Corporation 2024. Unauthorized use
prohibited. All rights reserved.

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© Neogen Corporation 2024. Toute utilisation non